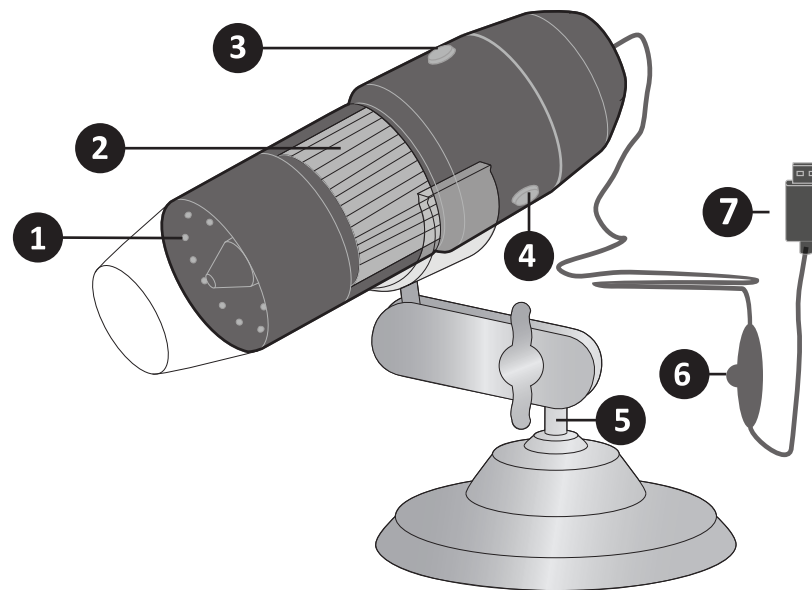
Microscope Numérique USB
USB Digital Microscope
USB Digitale Microscop
USB Digitalmikroskop

DESCRIPTION DU PRODUIT/PRODUCT DESCRIPTION/PRODUCT OMSCHRIJVING
/PRODUKTBESCHREIBUNG

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- 1. 8- LED light
 - 2. Focus wheel
 - 3. Magnification ratio (ZOOM)
 - 4. Snapshot/ Video
 - 5. Stand
 - 6. Led Lighting control
 - 7. USB line
- 1 CD ROMs included.

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SPECIFICATION

Image Sensor	1.3 Mega Pixels
Video capture resolution	Standard 640*480, Max 1600*1200
Focus Range	10mm ~ far view Manual focus from 10mm to 500mm
Magnification range	100X-200X, 50X-400X, 50X-500X, 50X-600X, 800X, 1000X
Video Format	AVI
Video encoding	MPEG2
Video Resolution	640*480mm or 1280*720mm
Video frame rate	30f/s under 600 LUX Brightness
Photo Format	JPEG
Picture resolution	160*120mm up to 8000*6000mm
Light Source	8 LED (adjustable by control wheel)
PC interface	USB 2.0
Power supply :	USB Port (5V DC)
Operation System	WINDOWS XP/VISTA/7/8 32/64 bit system
System Requirement:	Pentium Computer with 700M Hz & above, 20M HD Space CD ROM Driver, 64MB RAM, Direct X VGA Card

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DRIVER INSTALLATION PROCEDURE

1. Before connecting the microscope camera, turn on your PC and place the installation CD into the CD-ROM drive.

Note: Before proceeding with the installation, close the antivirus, in case it is opened, because the process of installation may be blocked by the antivirus software. In case it occurs, let the antivirus software allow all processes before installing.

2. Open "My computer" and double klik "the DVD/CD driver" to open the CD.

3. Click to install the driver and follow each step indicated below and make sure to click on each button circled in red in the following windows.

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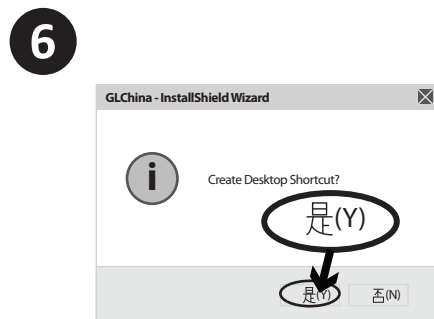
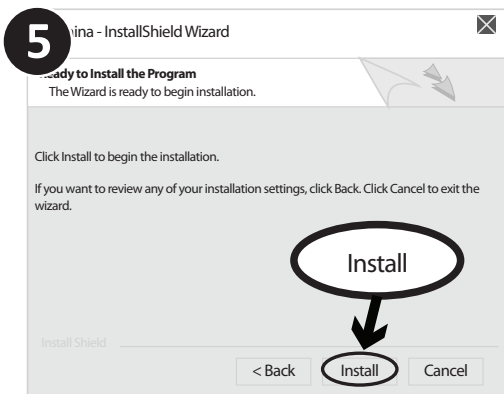
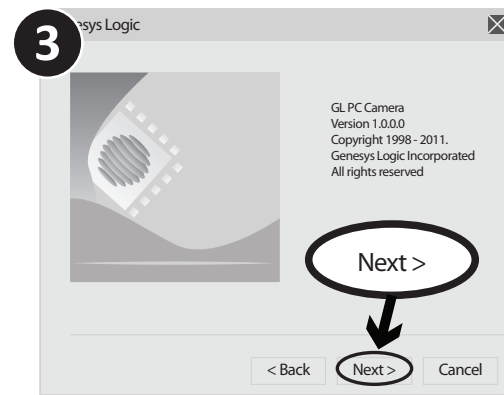
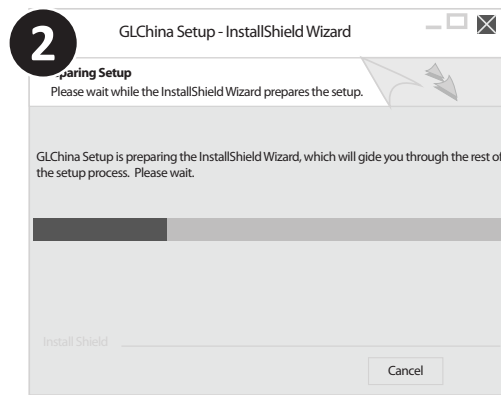
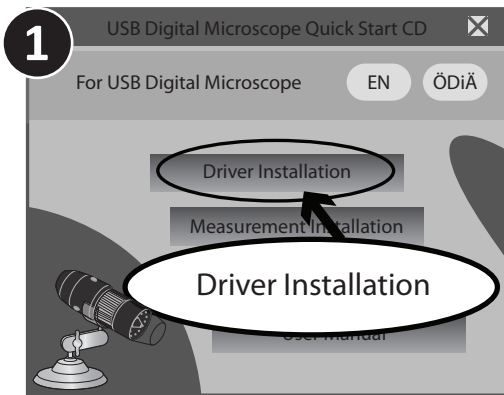
When installation is finished, reboot your computer at once if it is not rebooting. If the installation has succeeded, the following icon will be shown on the toolbar or the desktop.

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HOW TO USE YOUR MICROSCOPE CAMERA

1. Connect the USB cable of your microscope camera to the USB port of the computer.
2. Double click the camera device icon  on the desktop or toolbar. After clicking the interface below will show.

3. Hold your microscope camera in a vertical way to have a right image. Put it on to-be-observed object.
4. Adjust the brightness of the light with the LED lighting control (6). Then adjust the focus by rotating the focus wheel (2) with your thumb until the image is clear.
5. Take a picture by pressing either on the snap button or the icon .

Full Screen Display:

The real time display can be enlarged to full screen by clicking on the upper right corner of the window. To return from full screen mode, please click "Esc".

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
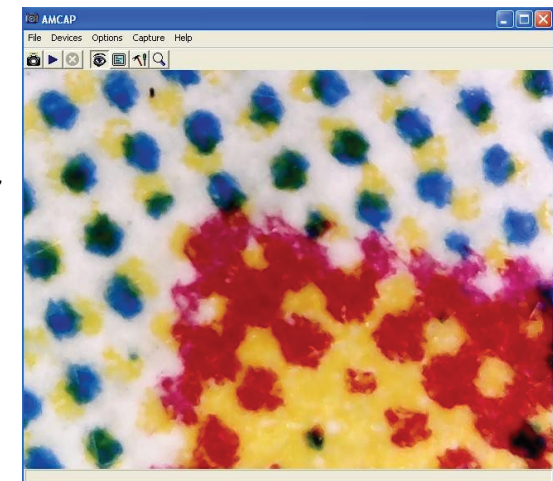
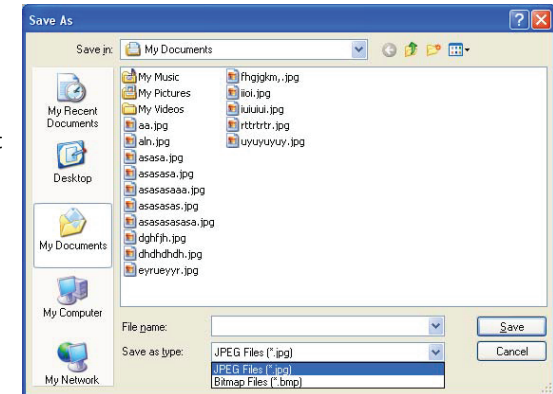
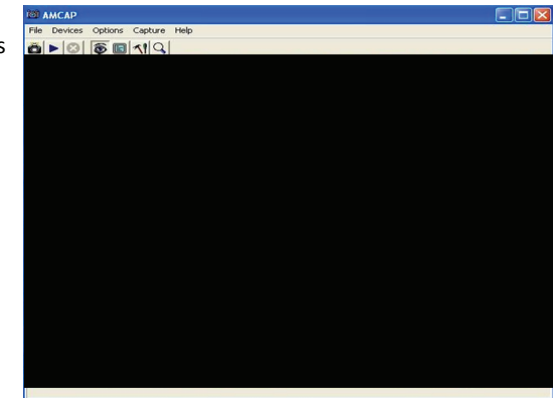
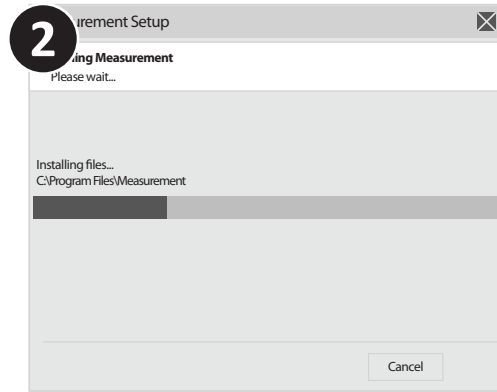
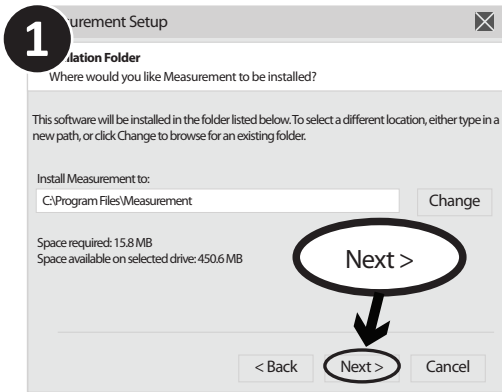
Kies en installeer de Meting

MEASUREMENT INSTALLATION

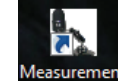
This function can only be installed on computers which operation system is Windows XP 32. Choose and install the Measurement

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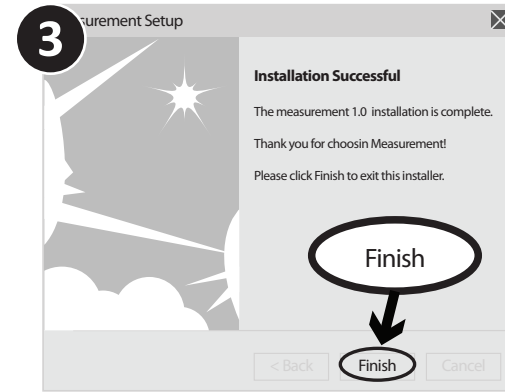Click "Finish" and the installation completes

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
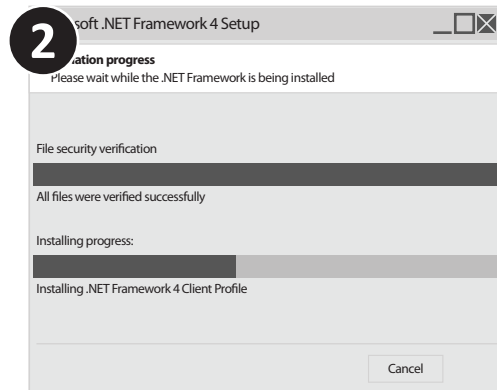
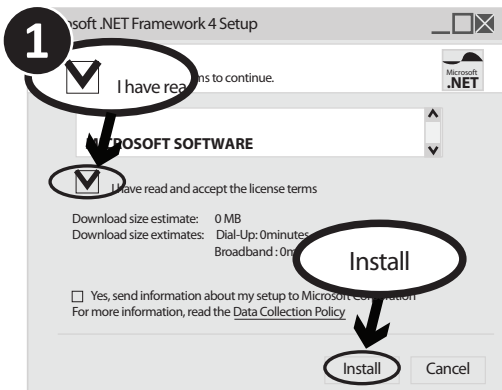
MICROSOFT .NET FRAMEWORK INSTALLATION

If the PC is Windows XP or older version, .Net Framework 4.0 is needed. Install it.

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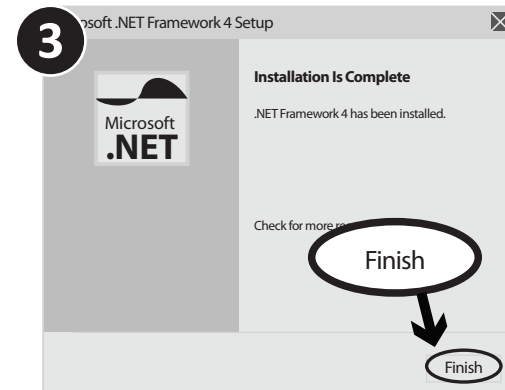FUNCTIONS OF OPTIONS

After the installation completes, double click at computer desktop to start using the software.

Note:

- Connect the camera to USB slot N°2 for using the video and photo capture.
- Remove the cap from camera to take the video and adjust the zoom to get a neat vision. Put back the cap to the camera when we use the microscope function.
- When we take the video, turn the camera's back upside for keeping the image straight.
- The software " measurement" works on the XP 32bit system. However, we can use the calibration plastic ruler (supplied) which has same functions as the software.

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Plug-in the devices, Display in operation window

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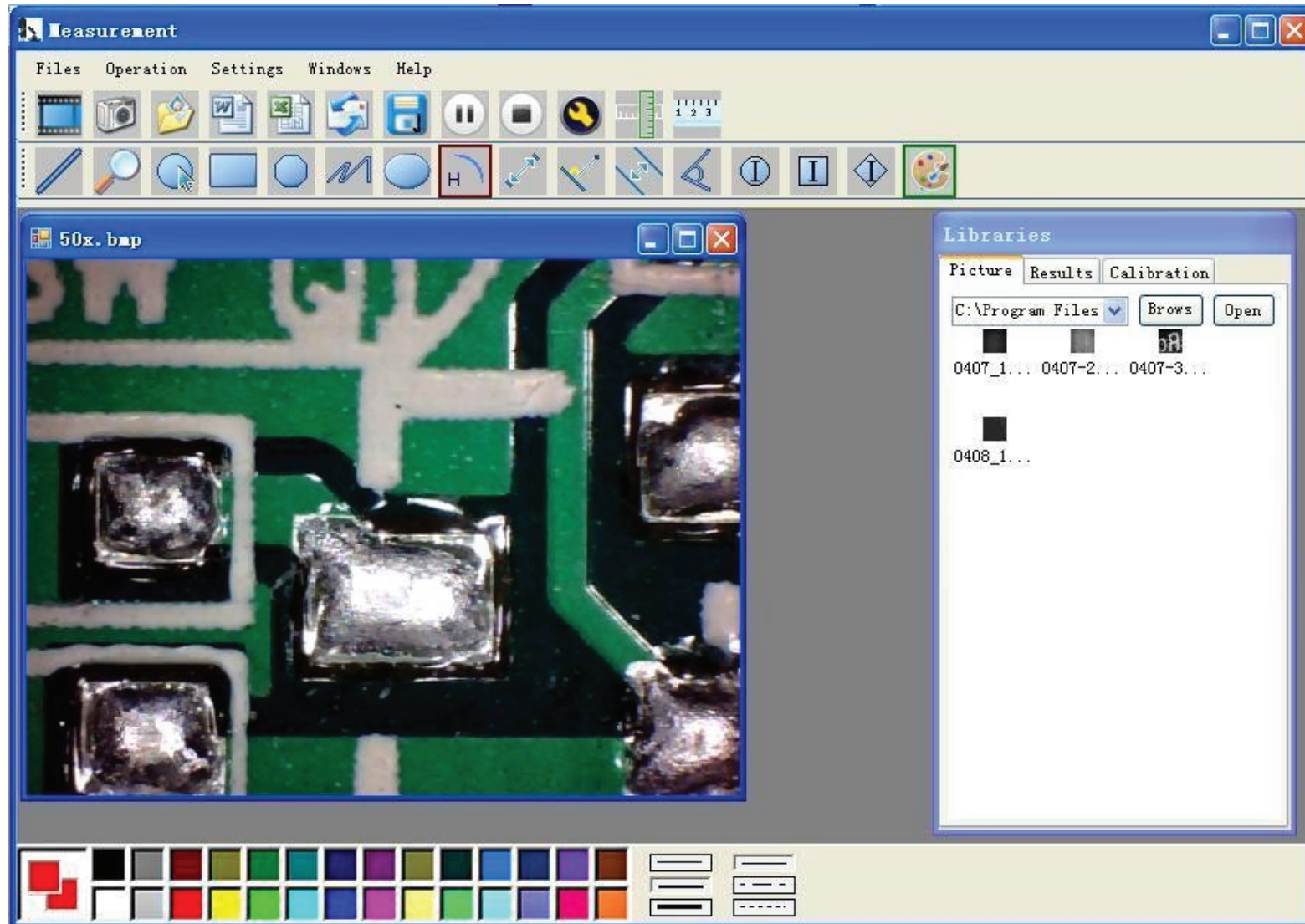
Choose the Image resolution from 320*240 to 1600*1200, default at VGA Mode640*480

Click  To use measurement function.

Click  to open pictures, and click  to save pictures

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

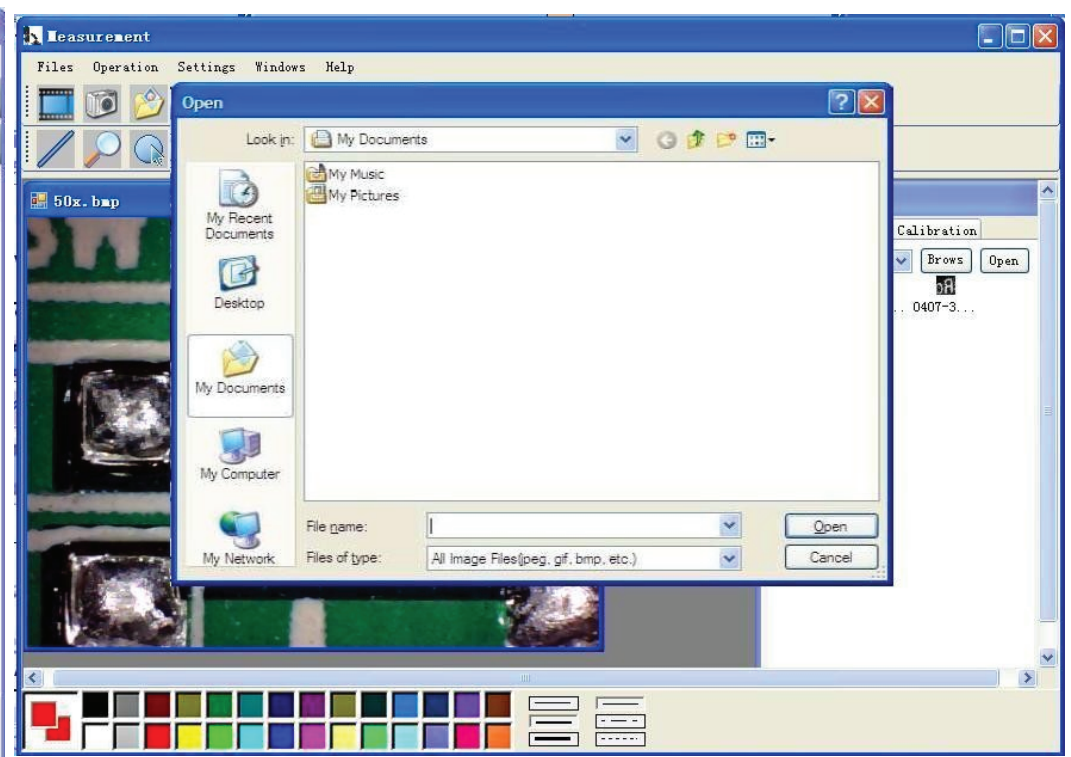
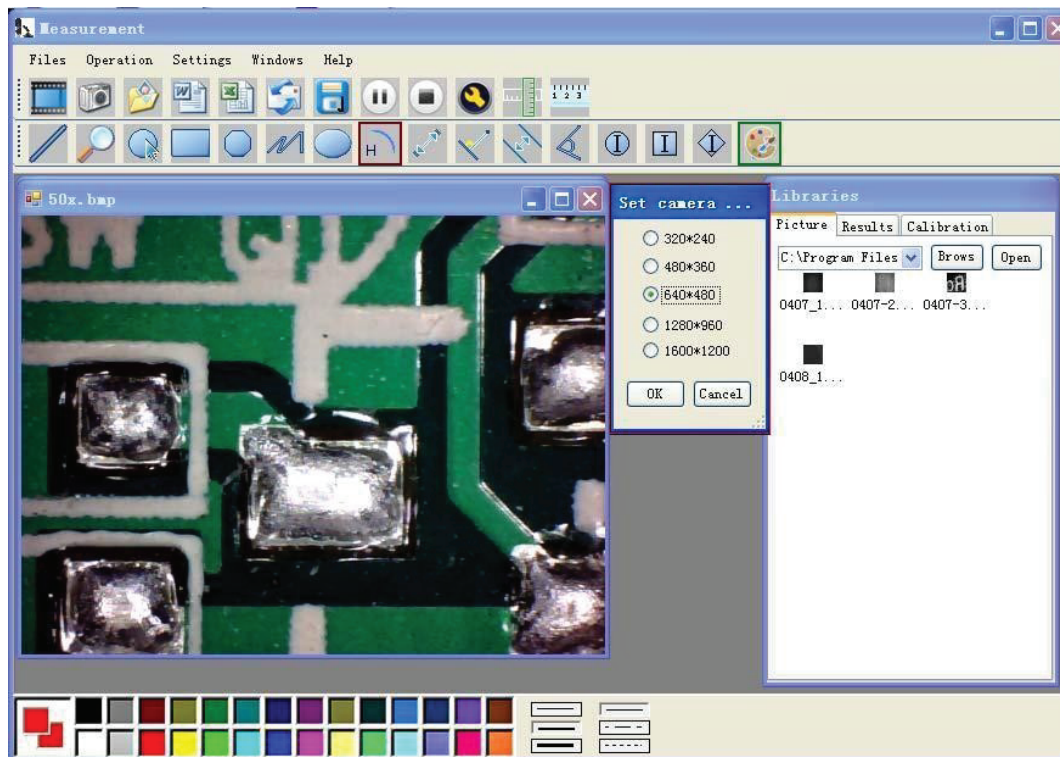
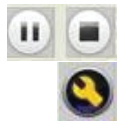
Vi deo PLAY (Pause), STOP

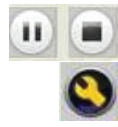Image Parameter including Image setting and special effects

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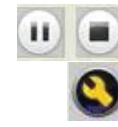Properties [Close]

Zoom | **Special Effect** | Video Proc Amp | Camera Control

Brightness		17	<input type="checkbox"/>
Contrast		100	<input type="checkbox"/>
Hue		0	<input type="checkbox"/>
Saturation		9	<input type="checkbox"/>
Sharpness		0	<input type="checkbox"/>
Gamma		22	<input type="checkbox"/>
White Balance			<input type="checkbox"/>
Backlight Comp		1	<input type="checkbox"/>
ColorEnable	<input type="checkbox"/>		

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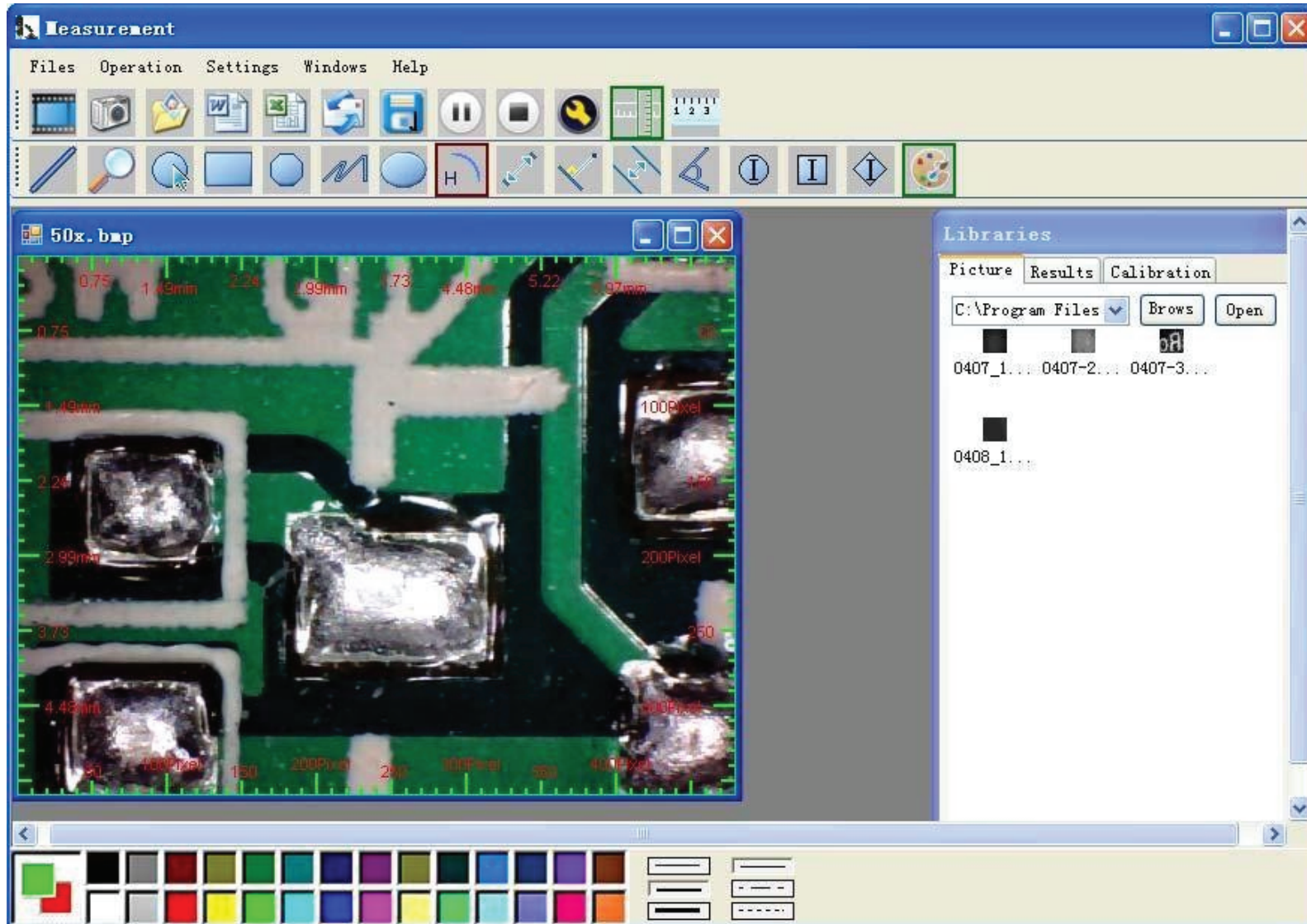
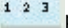
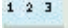
HOW TO DO THE CALIBRATION

Click  start calibration make sure certain accuracy. Our experience told us that measurement could be controlled at 99% high precision. Measure tolerance could be caused from human eye distortion and defocus. More practice is helpful to measurements.

Key in the Measure data (Pixel), Unit, and Click 'OK'.

Warning: Please do calibration again once Magnification Ratio changed.

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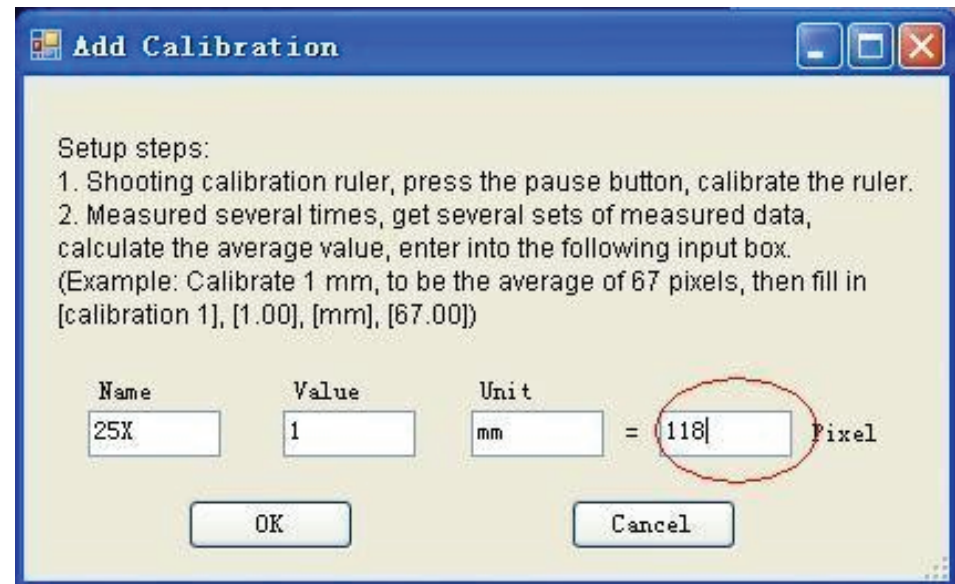
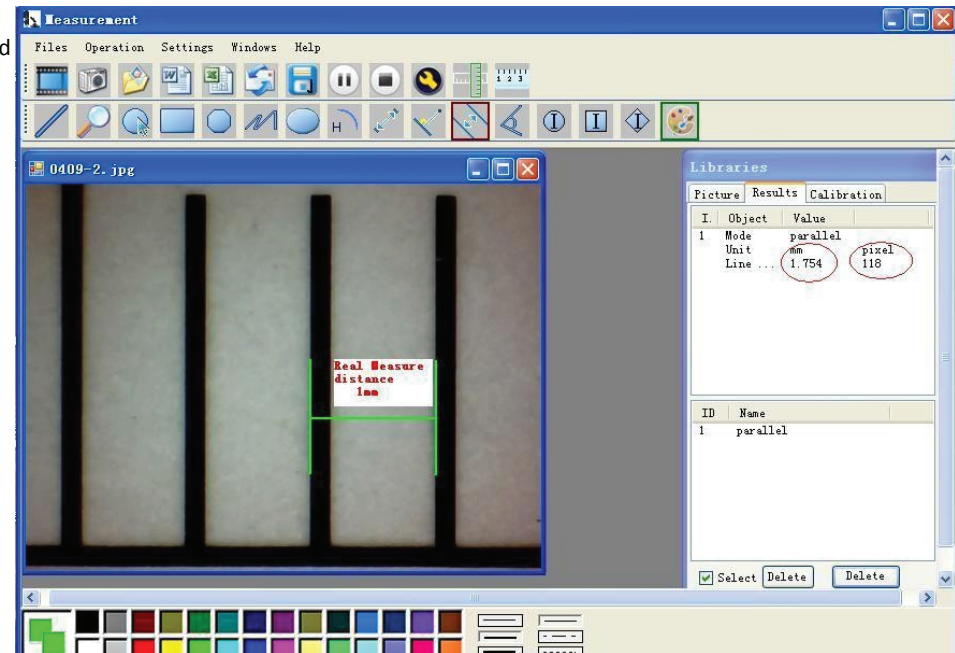
OTHER FUNCTIONS

Color setting of display character

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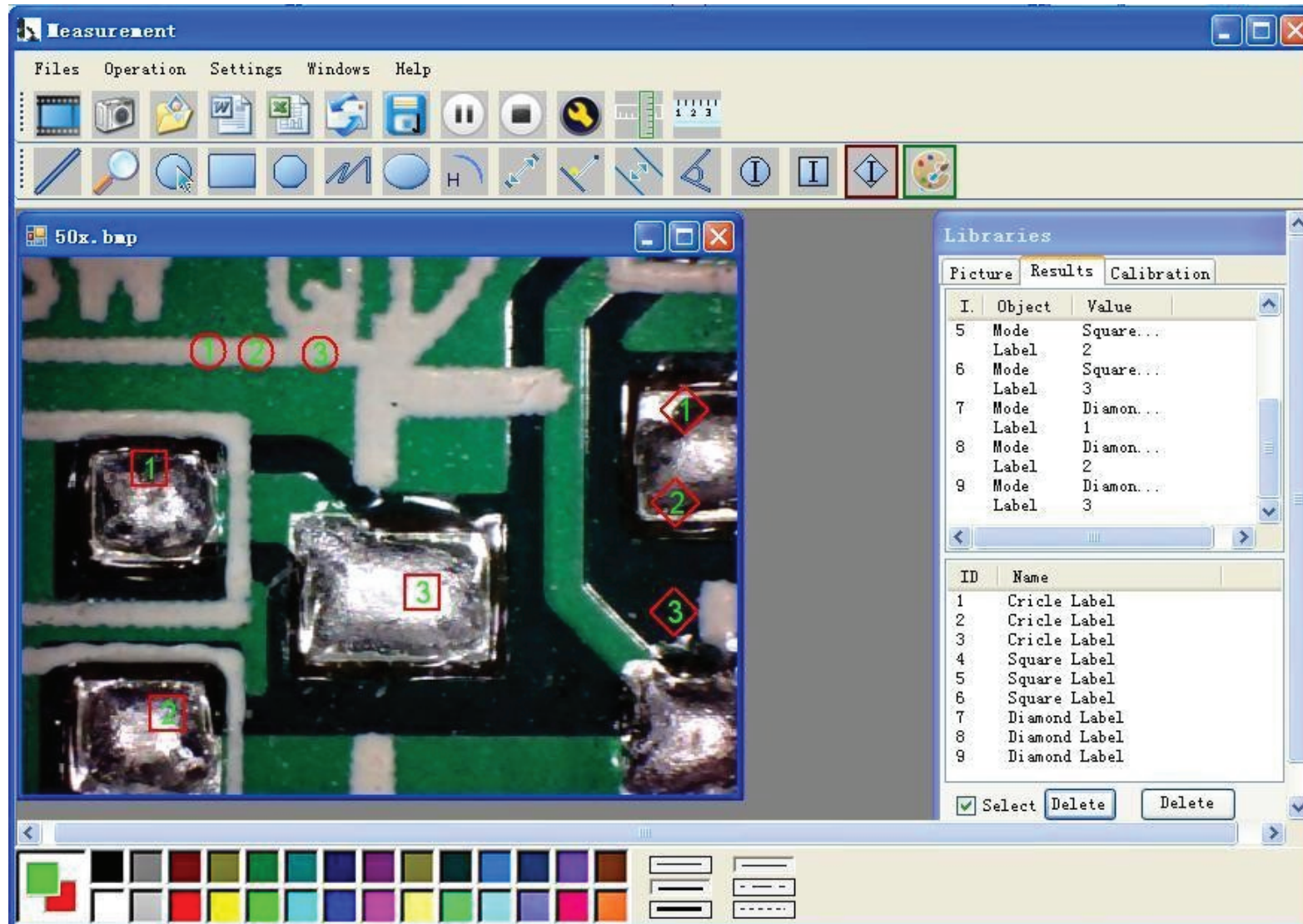
Example: Measure data under word file

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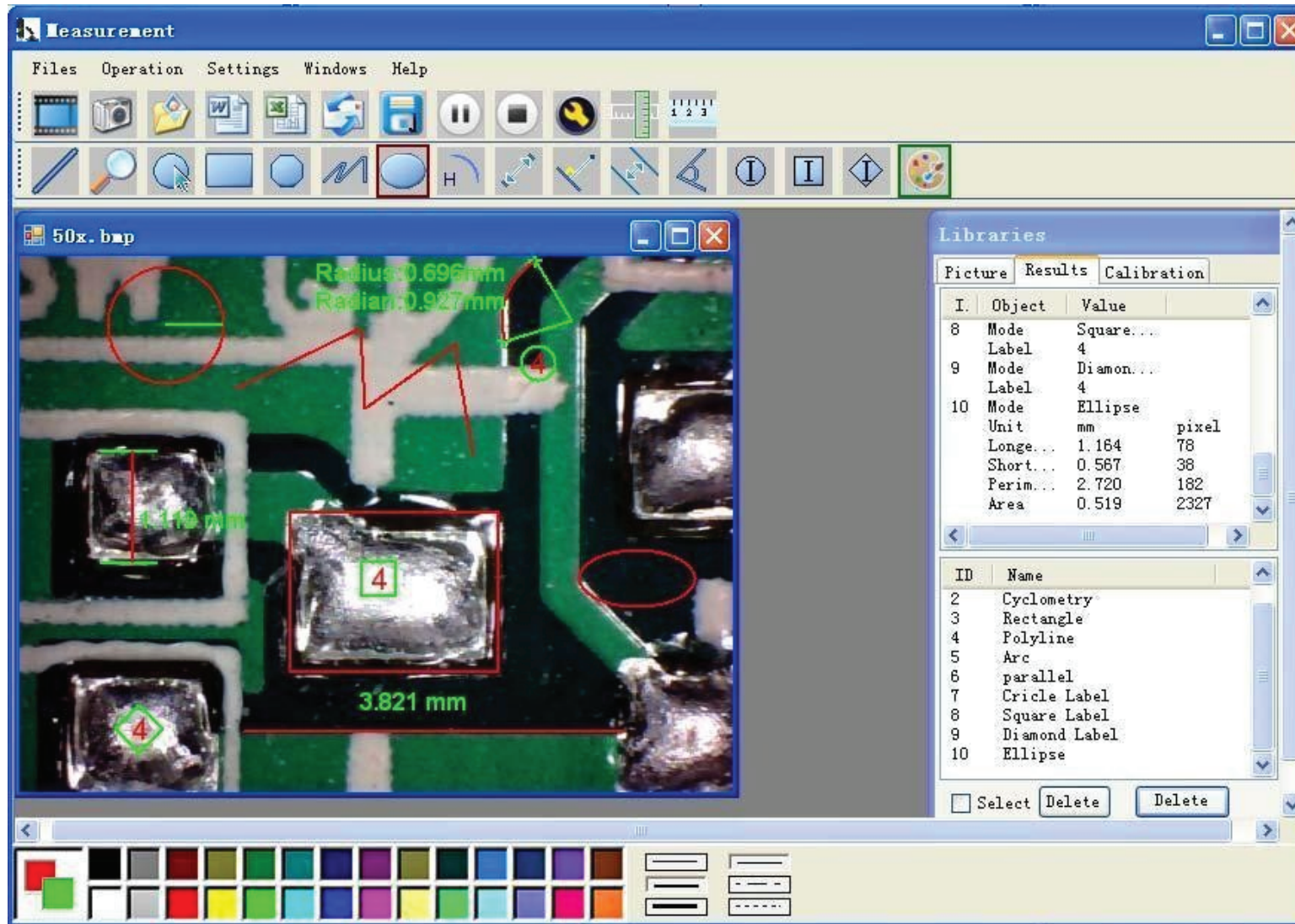
Line, ring, Square, Polygon, Set-line Length, Oval, Arc, Point to Point distance, Point to Line distance, Parallel line, Angle

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For example 1
Line measurement:

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

For example 2
Rectangular measurements:

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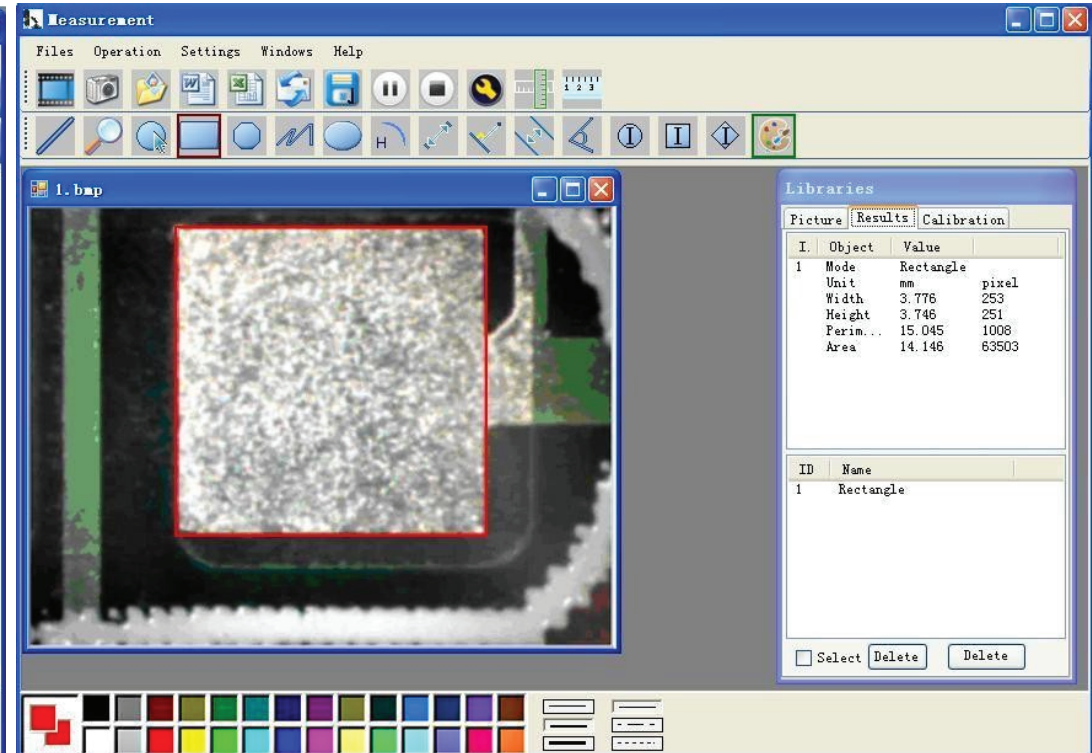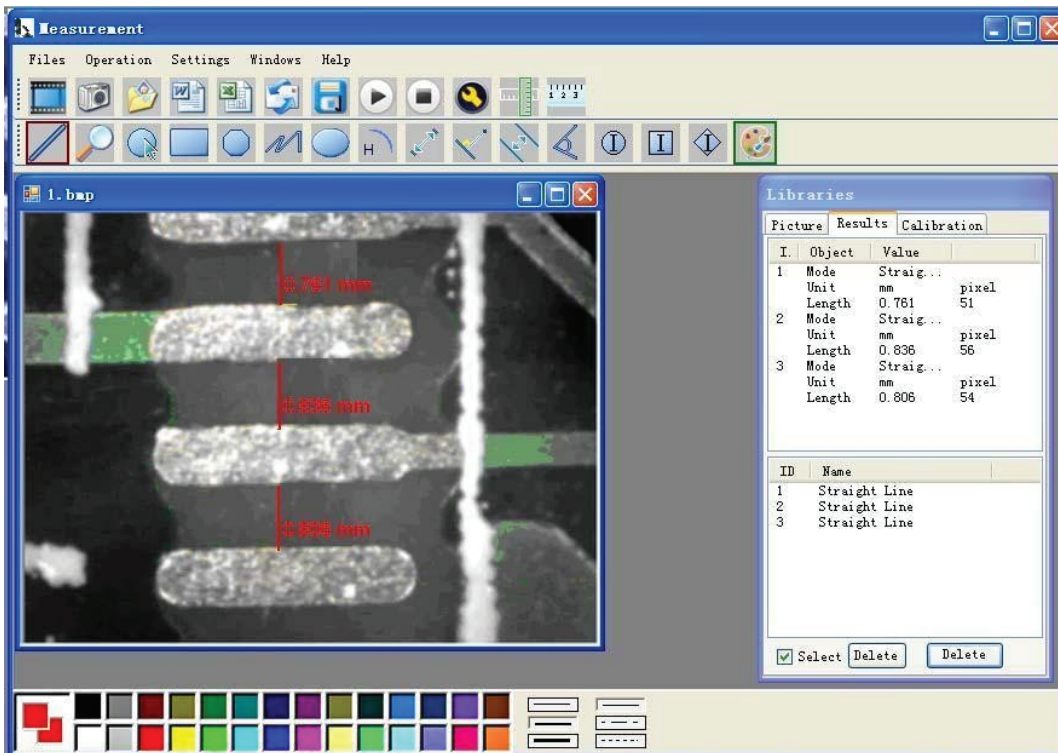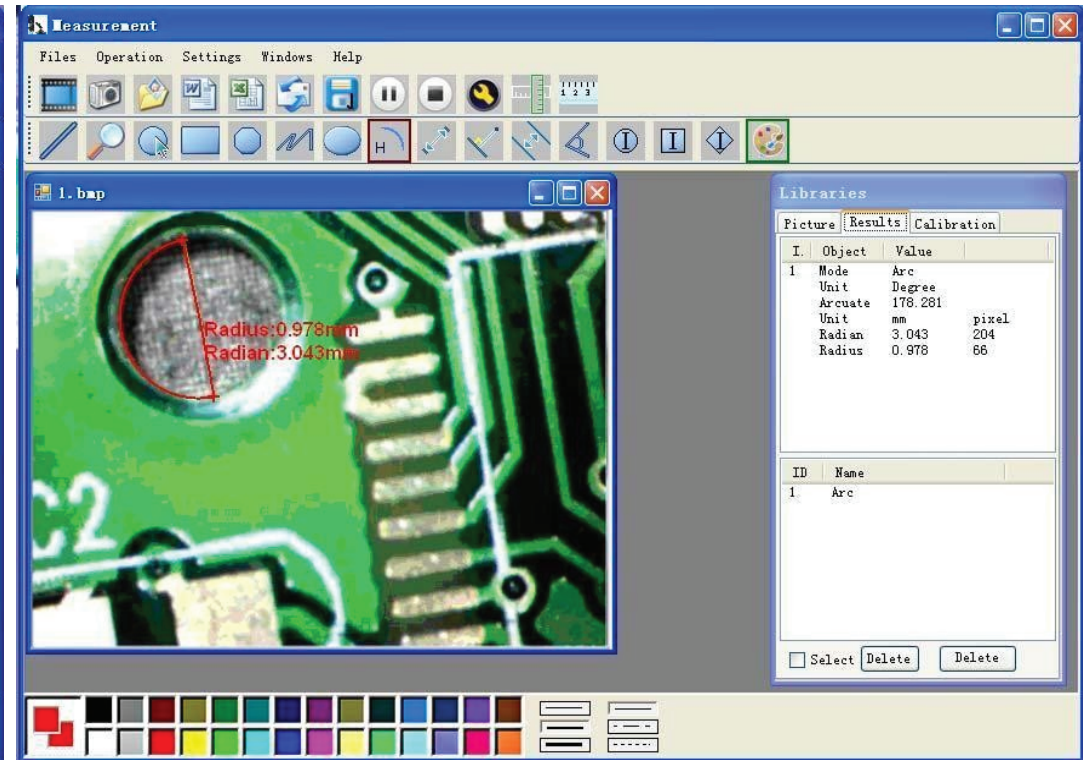
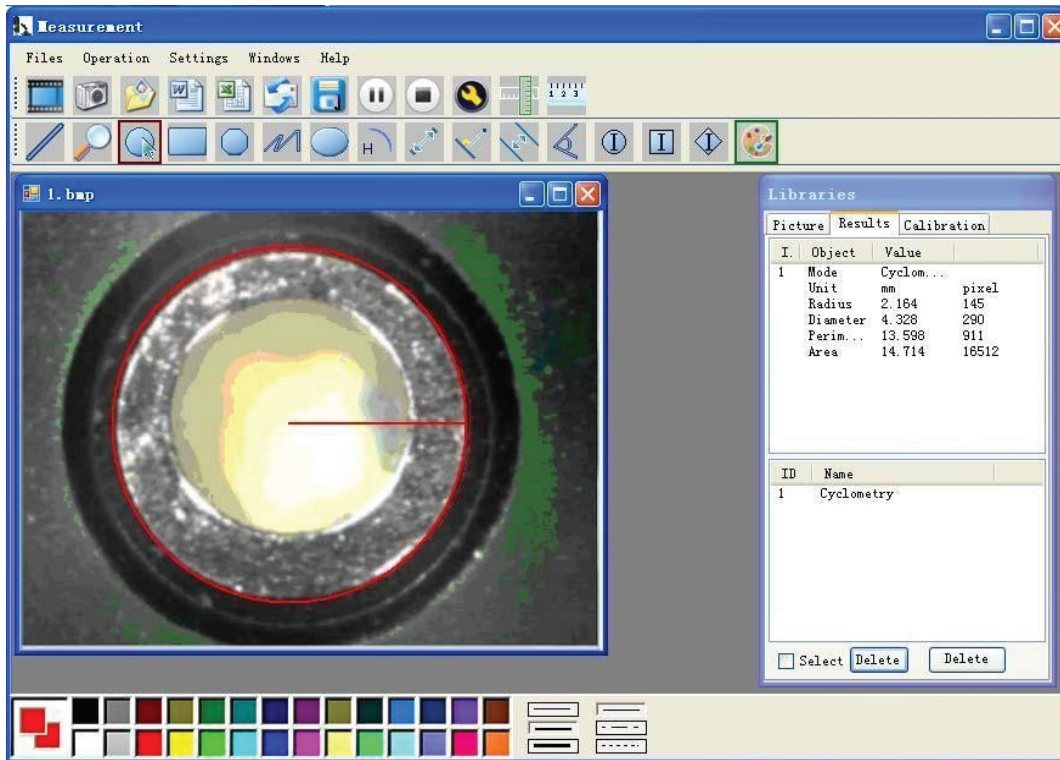
USAGE OF MICROSCOPE MICROMETER CALIBRATION RULER

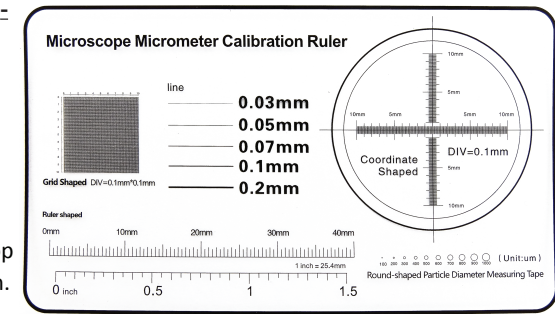
The microscope has slightly different magnification factors, the micrometer calibration ruler is used to calibrate your microscope for precise measurement. This ensures that you will be making accurate measurements with your microscope.

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WARNINGS

- o Do not shine light directly into eyes.
- o This is not a toy. Keep away from children.
- o Keep it away from heat sources, direct sunlight, humidity, water or other liquids.
- o Do not use it with wet hands.
- o Do not drop, short-circuit, modify or puncture it.
- o Do not disassemble the digital Microscope or change the interior parts, it can cause damage.
- o Do not touch the lens with your fingers.
- o In case the product gets wet, leave PC connection immediately.
- o Do not disassemble or dry by hair dryer. Send to repair center if the digit microscope was effected by liquid or other elements.
- o Do not allow this product to have contact with water or other liquids. If water or other liquids enter the product interior, immediately unplug the product from PC.
- o Continued use of the product may result in fire or an electrical shock.
- o Need to read User's Manual before using and storing it.
- o To avoid electric shock, remove all electricity source before moving PC.
- o Do not use any detergent or volatile solution, which will damage the camera case, painted surfaces & lens. Only use a soft cleaning cloth.

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WARNING: You should not dispose of this device with your household waste. A selective collection system for this type of product is implemented by your local authorities. Please contact your local authorities to find out how and where collection takes place. These restrictions apply because electrical and electronic devices contain dangerous substances that have harmful effects on the environment or on human health and must be recycled.

This symbol indicates that electrical and electronic devices are collected selectively. The symbol shows a waste container crossed out with an X symbol.

If the power cable is damaged, it must be replaced by the manufacturer, his after sales service, or a person qualified in this area in order to avoid any accidents arising.

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Importé par Prodis SAS, 1 rue de Rome, 93110 Rosny-sous-Bois, France/ Fabriqué en RPC
Imported by PRODIS SAS - 1, rue de Rome F-93110 Rosny-sous-Bois